

## FACT SHEET

# ● Bubblemeter



Bubblemeter 34 – 36 ml



Bubblemeter 500 – 560 ml

### **Available sizes:**

34 – 36 ml

40 – 80 ml

60 – 150 ml

500 - 560 ml

***others on request!***

In routine analysis the Burghart Bubblemeter is used to check the puff volume of smoking machines according to ISO 4387.

The operator sets the soap-film bubble to the zero position, by means of a pump and fine adjustment. Afterwards the connection tube of the Bubblemeter will be plugged into the smoke port of the smoking machine and a test puff is initiated. The soap-film bubble rises up with the air being drawn in. At the end position of the soap-film bubble the puff volume can be read off the graduation.

The Bubblemeter Burghart is easy and quickly to operate and offers high precision measuring results.

Supplied with a Factory Certificate. Optional and extra cost it will be supplied with a certificate for the volume graduation.

Subject to alterations.