

Bubblemeter



Bubblemeter 34 – 36 ml

Available sizes:

34 – 36 ml

40 – 80 ml

60 – 150 ml

others on request!

The Bubblemeter is a device to test the puff volume of a smoking machine in accordance with ISO 4387.

The wetting agent solution must be filled into the Bubblemeter to generate a soap-film bubble in the burette.

The operator sets the soap-film bubble, by means of an inflation bulb, to the zero position. Afterwards the connection tube of the Bubblemeter will be plugged into the smoke port of the smoking machine and a test puff is initiated. The soap-film bubble rises up until the test puff is finished. The distance between the zero position and the current position of the soap-film bubble is the puff volume.

Subject to alterations.