

Bubblemeter



Bubblemeter 34 – 36 ml

Available sizes:

34 – 36 ml

40 – 80 ml

60 – 150 ml

others on request!

The Bubblemeter is a device to test the puff volume of a smoking machine in accordance with ISO 4387.

The wetting agent solution must be filled into the Bubblemeter to generate a soap-film bubble in the burette.

The operator sets the soap-film bubble, by means of the syringe, to the zero position. Afterwards the connection tube of the Bubblemeter will be plugged into the smoke port of the smoking machine and a test puff is initiated. The soap-film bubble rises up until the test puff is finished. The distance between the zero position and the current position of the soap-film bubble is the puff volume.

Subject to alterations.